


# PASPort™

## USB Link

PS-2100A

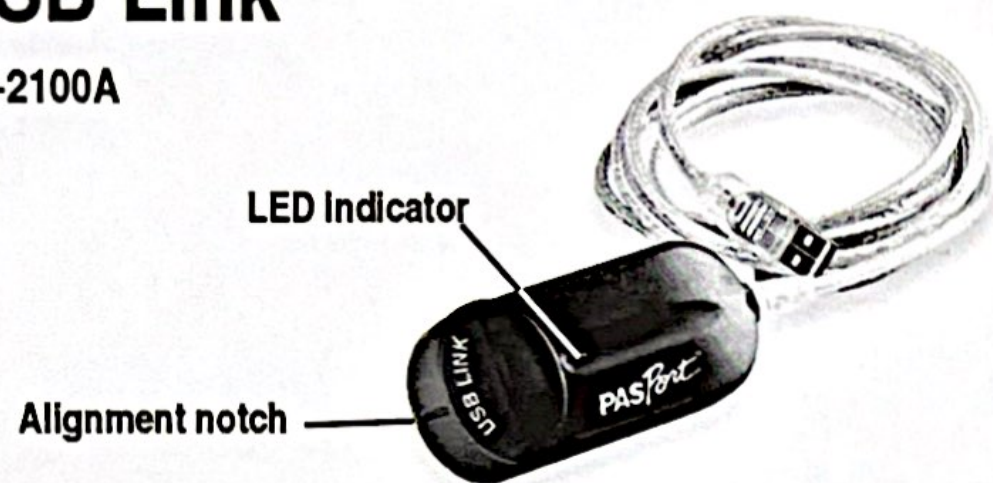

### USB Link Specifications

<b>LED ON:</b>	Indicates the USB Link is connected and software recognizes the connection.
<b>LED OFF:</b>	Reconnect the USB Link to the USB port or hub.
<b>One-way Connection:</b>	Align the notch on the USB Link with the groove on the sensor or sensor plug.
<b>Hot Plugging:</b>	USB Links can be connected to or disconnected from a USB port without turning off the computer.

## USB Link Quick Start

The PS-2100A USB Link is a simple PASPORT Link Device that allows PASPORT sensors to communicate with a USB-compatible computer.

The Universal Serial Bus (USB) is a computer protocol that allows advanced communications between peripherals (e.g. sensors, keyboards, mouse devices, monitors, printers, etc.) and your computer.

### Computer Requirements

- At least one USB port
- EZscreen or DataStudio Software (version 1.5 or later)
- **Macintosh** - Mac OS 8.6 or higher
- **Windows** - Windows 98 or 2000

### Equipment Setup

1. Install the EZscreen software included with all PASPORT sensors. EZscreen software only needs to be installed once.
2. Connect a PASPORT sensor or sensor plug to the USB Link.
3. Connect the USB Link to a USB port on your computer or USB hub. (Look for the USB icon shown to the right.)
4. The software launches when it detects a PASPORT sensor. Select **EZscreen** or **DataStudio**.

